


Instructions for Whirl-Pak® THIO-BAGS®

Caution: Do not allow the thio tablet to fall out of the bag.


OPENING



Tear off top along perforation.

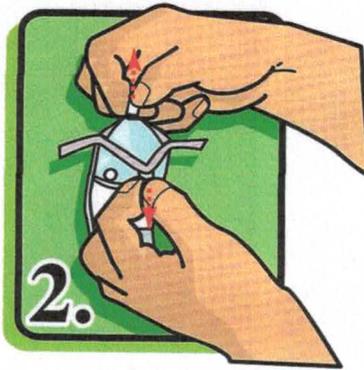

Pull tabs outward to open bag. Sometimes a pull on the bottom is also helpful.

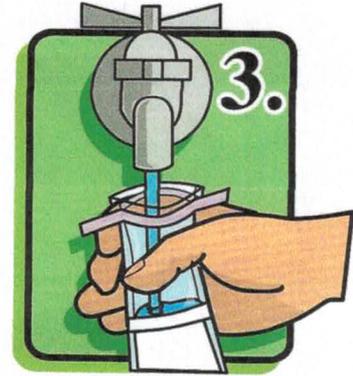

Fill bag to required fill line. The 4 oz. fill line will ensure that 100 mls are collected.

CLOSING



Pull wire ends to close bag.



Whirl bag three complete revolutions. (Do not roll tapes down.) Leave airspace for mixing in lab.



Turn tape inward on opposite face of fold.

NASCO • FORT ATKINSON, WISCONSIN • MODESTO, CALIFORNIA SPECTRUM/NASCO • NEWMARKET, ONTARIO, CANADA
PRINTED IN U.S.A. U05352 /RV 9-09

Instructions for Dropping off Well Water Sample:

WITH A BALL POINT PEN, WRITE YOUR NAME, DATE AND LOCATION YOU TOOK THE SAMPLE (OUTSIDE FAUCET, ETC.)

DO NOT TOUCH TOP OF BAG TO FAUCET OR WITH FINGERS AFTER TEARING OFF PERFORATION.

PUT THE SAMPLE IN A COOLER WITH AN ICE PAK AND TAKE TO: THE WATER SPIGOT, 5806 E HIGHWAY 22, PANAMA CITY FL. PHONE: (850) 871-1900

THE SAMPLE SHOULD BE SUBMITTED WITHIN FOUR HOURS OF BEING TAKEN.